WiseCircu[®] WCB

Umwälzthermostat

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Circulation Water Bath

Operating Manual

gilt für / written for
WCB -6/ -11/ -22

witeg Labortechnik GmbH

(English manual at page 13)

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Contents

1 Introduction	14
2 Safety Instructions	14
2.1 Description of symbols.....	14
2.2 Instructions	14
3 Product description	16
3.1 Key features	16
3.2 General survey.....	17
3.3 Operating surface.....	18
3.4 Term description.....	18
4 Packing content	19
4.1 Devices with bath	19
4.2 Immersion devies	19
5 Positioning the device and commissioning	19
6 Job description	20
6.1 Menu structure.....	20
6.3 Set turn-on delay	20
6.4 Set temperature	21
6.5 Justification	21
6.6 Key lock.....	21
6.7 Storage function	21
7 Maintenance and cleaning	22
8 Troubleshooting Guide	22
9 Technical data	23
10 Certificate	24

1 Introduction

Thank you for purchasing our **WiseCircu® WCB Bath Circulator**. You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- It is important that the device stands up on a horizontal, stabile and firm subsurface to avoid vibrations.
- If you set the device in a site with vibrations, make sure that the device can not slip away.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in area with potentially explosive atmosphere.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.
- Make sure, that the safety switch is adjusted up to 30°C over max. temperature. Otherwise the heating maybe doesn't work.

- Do not put heat-sensitive items at the device.
- Warn all attendees if you leave the device unattended.
- Pay attention when you lift up the lid, hot condensation water may drop down or hot steam could come out.

- Do not clean the scorching device.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

3 Product description

3.1 Key features



The **Wisdom WiseCircu® WCB Circulation Water Bath** is designed for chemical, biological, medical and pharmaceutical applications. The device was developed for simplest usage and a very long life.

Key features are as follows:


- Simplest operating by MODE button and Jog-Shuttle knob.
- Fast commissioning.
- Long life by simply and solid construction.
- Highest temperature accuracy by digital fuzzy controller.
- Temperature and delay time showed on illuminated LCD.
- Storage function for set time and temperature values.
- Turn-off an -on delay from one minute up to over 99 hours in one minute steps adjustable.
- Signal sound after end of turn-off delay.
- Easy to justify.
- Key lock against inadvertent changing of set speed and timer values.
- Over temperature protection by safety switch.
- A lot of laboratory devices are mountable.

3.2 General survey

The functionality of the described components is the same at each type of the **WCB**-series. Only the position may differ from type to type.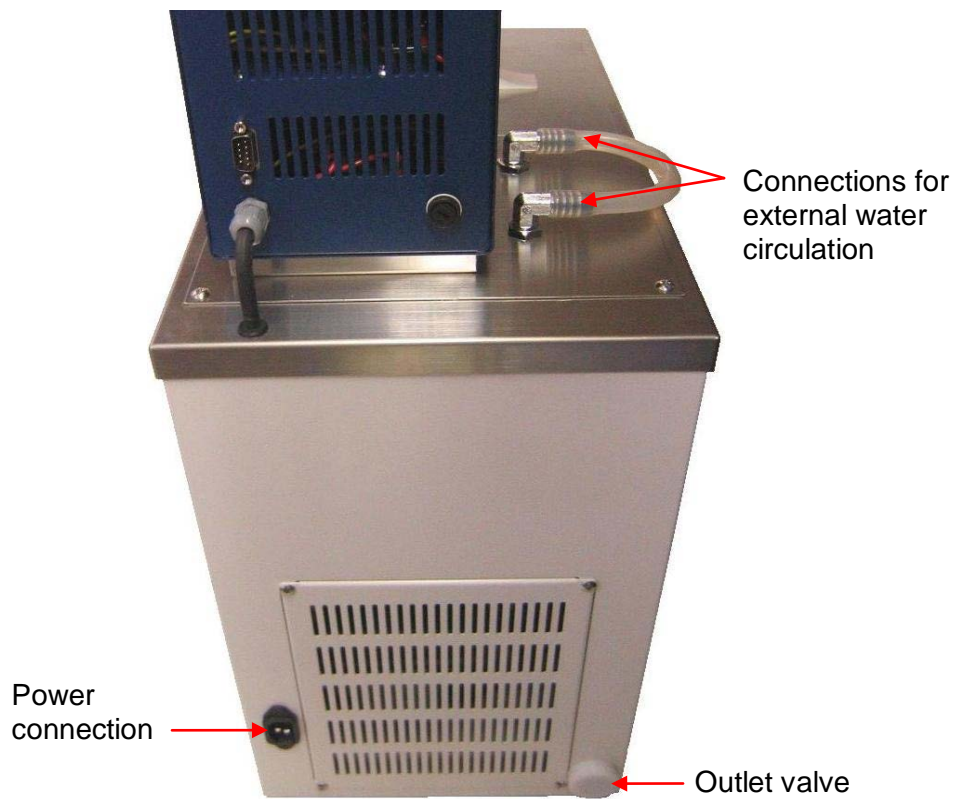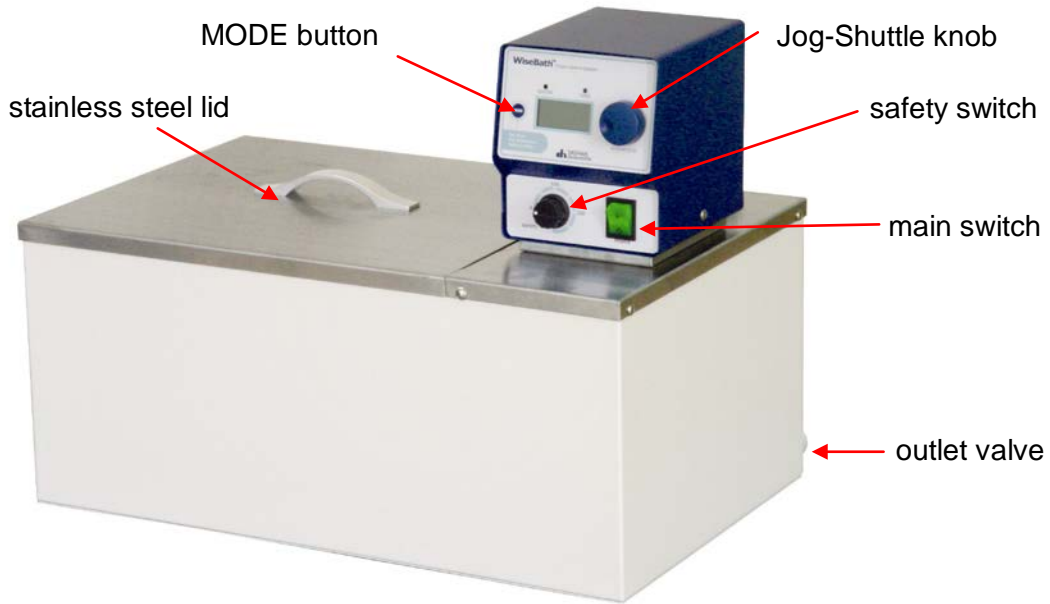

3.3 Operating surface

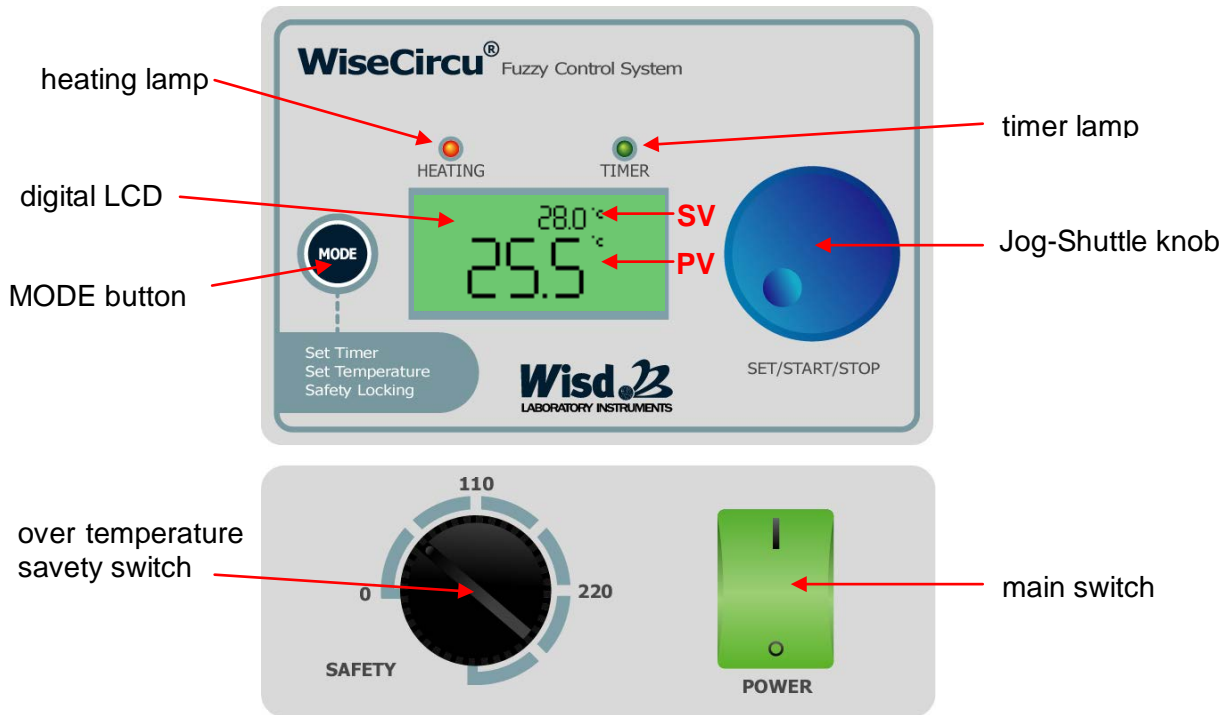3.4 Term description

SV: set value (upper line, in key lock menu no SV)

PV: present value, measured by the device (lower line, in key lock menu also upper line)

Main switch

- On / off switch.

Timer lamp

- Is blinking in menu items for timer modulation
- Shines continuous in the key lock menu item.

Heating lamp

- Shines if the heater is activated.

Jog-Shuttle knob

- Changes values by turning and stores, confirms, stops and starts by pushing.

Safety switch

- The safety switch turns the heating off if your set maximal temperature is overshoot.
- Set the maximal temperature roughly 10°C-20°C over the desired maximal temperature, otherwise the device maybe not reaches your set value.

MODE button

- For changing between the different menu items and the key combination.

Key combination

- To reach the sub menu Items you first have to hold the Mode button pushed and push in addition the Jog-Shuttle knob.

Connections for external water circulation

- To reach the sub menu Items you first have to hold the Mode button pushed and push in addition the Jog-Shuttle knob.

4 Packing content

4.1 Devices with bath

quantity	DH.WCB00106	DH.WCB00111	DH.WCB00122
1	WCB-6	WCB-11	WCB-22
1	connection cable		
1	stainless steel lid		
1	tube (ca. 10 cm)		
1	operating manual		

4.2 Immersion devies

quantity	DH.WCB10011	DH.WCB10022
1	WCB-11H	WCB-22H
1	connection cable	
1	operating manual	


The immersion devices does have the same performance features like the bath-devices **WCB-11** and **WCB-22**.

5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol: 
2. Remove the package and perhaps soiling carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your costumer service.
4. Set the protection switch 30 °C over desired bath temperature otherwise the heating coil does not work.
5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
6. Make sure that you follow all safety instructions.
7. Connect the device with the power outlet.
8. Fill the operating liquid in.
9. Switch the device on and start the process (look at job description).

6 Job description

6.1 Menu structure

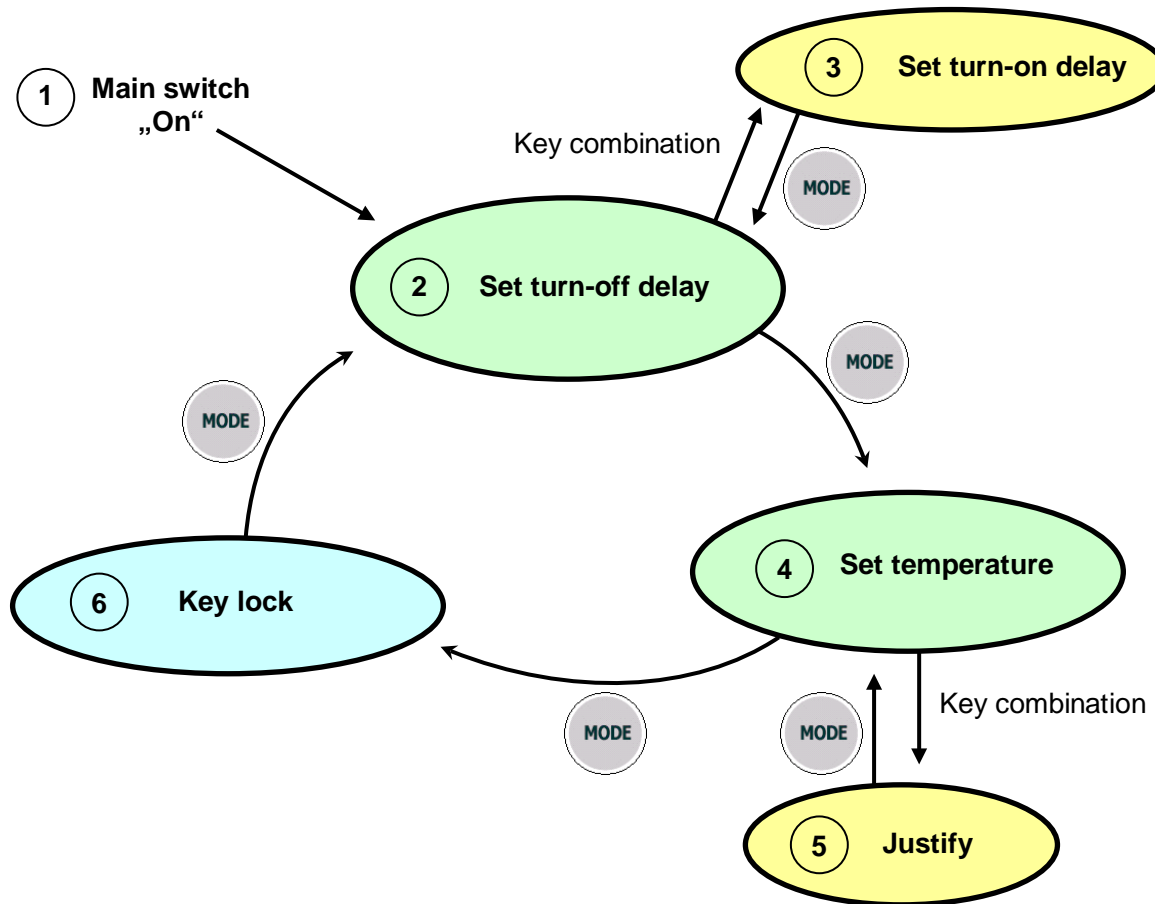6.2 Set turn-off delay

The turn-off delay timer serves as automatically stop function. It turns off the heating after the timeframe you set is elapsed.

- Use Mode button to reach next menu without setting a delay time.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- Turn the Jog-Shuttle knob to set your delay time. Push the Jog-Shuttle if you reached the desired value to confirm it. With the confirmation the timeframe starts elapsing.
- If you want to change the value after confirmation, you have to push the Jog-Shuttle once again. The PV initiates blinking and you can modify the value again.
- To confirm the value and start the delay time push the Jog-Shuttle yet again.

6.3 Set turn-on delay

The turn-on delay timer serves as automatically start function. It turns on the heating after elapsing of the timeframe you set.

- To reach this sub menu item you have use the key combination (look at 3.3) in the menu “set turn-off delay”. **LCD shows “d-tr”**.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- Turn the Jog-Shuttle to choose your desired value and probate it by pushing the

Jog-Shuttle. The delay time starts to run by probating the value.

- If you want to change the confirmed value, you have to push the Jog-Shuttle once again.
- Turn the Shuttle knob clockwise to increase the time and counter-clockwise to reduce it.
- The turn-off delay time will not elapse before the turn-on delay time is executed.

6.4 Set temperature

- When you push the MODE button in the menu item “set turn-off delay” you reach the menu item “set temperature”. You can recognize this menu at “°C” in SV and PV line.
- The temperature is adjustable from 15 °C till 100 °C in 0,1 °C-steps.
- Push the MODE button once again if you want to work without the heating function. Then you reach the next menu (look at 6.1) without collimated temperature value.
- Choose your desired temperature by turning the Jog-Shuttle and confirm the set value by pushing the Jog-Shuttle.
- For changing the confirmed value again, you first have to push the Jog-Shuttle knob again. “HEATING” starts blinking at the LCD and the temperature is adjustable again.
- When the heating lamp is shining, the heater is activated.

6.5 Justification

If the temperature measured by you differs from the temperature the device is showing (PV) you have the possibility to compensate the divergence by the Offset. So you can balance differences of about ± 10 °C.

The Offset is calculated as follows:

Temperature measured by you – showed temperature (PV) = Offset

- To reach this menu you have to use the key combination in the menu item “set temperature”.
- Set the calculated value by turning the Jog-Shuttle. The Offset will be added to the PV immediately.
- Save the offset by holding the Jog-Shuttle pushed until the values in PV and SV line starts blinking.
- Push the MODE button to leave the justification menu.

6.6 Key lock

- To avoid accidental changing of the set values, push the MODE button in the menu item “set temperature”.
- In this menu item can you see the actually temperature in the SV-line and the remaining time of turn-off delay in the PV-line. The function of the Jog-Shuttle knob is deactivated.
- Press the MODE button once more to leave the key lock. Than you are in the menu item “set turn-off delay”.

6.7 Storage function

The **WiseCircu® WCB Circulation Water Bath** includes a storage function for temperature and time values. That is practical if you want to work with the same parameters often. Then you did not need to set them in every usage.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.
- Any stored parameters are erasable. Turn the value with the Jug-Shuttle till “0” and then hold the knob pressed for two seconds.

- If you want to overwrite a value it is not necessary to erase the old one. Set the desired value and use the Jog-Shuttle knob for storing.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> ✓ Check the supplied voltage is present. ✓ Check the power connection and the power cord for proper connection and integrity. ✓ Check if the MAIN Power Switch is "ON". ✓ Fuse is blown.
Unit Not Heating	<ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check if the set temperature (SV) confirmed with the Jog-Shuttle knob (by turn the Jog-Shuttle knob cannot change the set temperature – see point 6.4) ✓ Check the Safety switch is set higher (min. 10°C-20°C) than the set temperature (SV).
Heating, but unit not reaching temperature	<ul style="list-style-type: none"> ✓ Check the Safety switch is set higher (min. 10°C-20°C) than the set temperature (SV). ✓ Problem in the sensor or main controller.
ERROR 1	<ul style="list-style-type: none"> ✓ Temperature Sensor Error.
ERROR 2	<ul style="list-style-type: none"> ✓ Low level error. Please fill water in.

If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.

Reparations and amendments should only made by [witeg](#) authorized people. Contraventions exclude the warranty.

9 Technical data

		WCB-6	WCB-11	WCB-22
Dimension (BxTxH)	Bath (mm)	300 x 155 x 150	302 x 240 x 150	500 x 295 x 150
	External (mm)	330 x 185 x 385	338 x 275 x 385	540 x 335 x 410
Capacity		6 Lit.	11 Lit.	22 Lit.
Heating power		800 W	1200 W	2000 W
Temp.	Range	Ambient + 5 °C - 100 °C		
	Accuracy	± 0,1 °C		
	Uniformity	± 0,2 °C		
	Sensor	PT100		
Pump		14 Lit./min		
Controller		Digital fuzzy controller with Jog-Shuttle knob (Turn + Push)		
Display		Illuminated LCD		
Timer		99 h 59 min		
Safety device		Over temp. and over current protection, low filling level detector		
Material	Internal	Stainless steel		
	Capsule	Powder coated steel		
	Lid	Stainless steel		
Others		<ul style="list-style-type: none"> - Storage function (temperature and timer), - Key lock (Jog-Shuttle input disabled) - Alarm (sensor error, low filling level and timer-end) 		
Power supply		AC 230 V, 1N~, 50/60 Hz		

10 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	WiseCircu® WCB Circulation Water Bath
Typ	
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de

Version 1.6.25112013