# MINISPENSOR

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## Operation Manual MINISPENSOR

Please follow the step by step guide.

### General safety precautions:

Highest safety precautions should be used when dispensing corrosive, poisonous, radioactive or hazardous chemicals.

- Observe general safety regulations (e.g. wear safety goggles, protective clothing, etc).
- Carefully follow the operation manual.
- Use the instrument only with regard to resistance of material and for the purpose for which it is intended. In case of doubt concerning the suitability please contact the manufacturer.
- Always check the dispenser for leaks. Make sure that the suction and ejection canula are firmly in position before starting to use the instrument.
- NEVER use force! Breakage of any part will lead to hazardous exposition to the user/ operator as well as others.
- We only consider ourselves responsible in regard to safety, reliability and performance of the MINISPENSOR if:
  - ◆ The dispenser has been used in accordance to the instruction manual
  - ◆ Only the parts supplied by the manufacturer have been used

### Before operating the dispenser ...

Check if dispenser has not been damaged during transport.

The Minispensor is suitable for practically all liquids and dispensing methods.

#### EXCEPTION:

Solutions of hydrofluoric acid or of substances which form crystals, contain or form solids, as well as solutions of substances which are catalyzed by platinum and iridium alloy. Inorganic oxidizable solutions (e.g. biuret reagent) may precipitate metallic oxides, therefore clean the dispenser daily:

The temperature of the instrument and the reagent should not exceed 40 °C (104 °F).

The user has to ensure the compatibility of the dispenser and its chemical resistance for himself.

Select the appropriate length of the canula according to the bottle which will be used and push it on tightly. Screw the ejection canula on firmly.

### Do not use the plunger before the dispenser has been assembled completely!

The Minispensor fits to the corresponding 100 ml bottle, also supplied with an adapter A28/32; more adapters, see SPARE PARTS LIST

**Warning:** Avoid personal injury from chemicals. To protect yourself always wear eye protection and other appropriate safety equipment and clothing. Please follow all safety instructions and observe the operating procedures.

Safe and proper operation and function is only possible with the ejection canula as originally supplied. Never use damaged or deformed canulas. Do not use any other type of tubing. Regularly inspect the tubes for leakage and replace immediately damaged or deformed ones.

### Dispensing

Which fluid do you wish to dispense?

Be sure to observe all safety instructions and operating limitations.

Place the collecting vessel under the ejection tube. The ejection canula must face away from operator or any other persons at all times.

**Warning:** For your own personal safety:

Never press down the plunger as long as there is no vessel to collect the liquid.

**CAUTION:** If the valve is damaged, reagent may drop out of the ejection canula!

### Function and ventilation

Different than other dispensers Minispensor is the only dispenser which enables a non-tiredsome use. Long series of work can be easily done. This new innovated dispenser can be operated with only three fingers. The plunger has to be pushed down with the thumb (just like with an automatic pipette), the selected volume will then be obtained through a spring.

Before the first use the dispenser has to be ventilated. The air caught in the supply pumps can be removed with 2-3 pumps

### Volume adjustment

The Minispensor is supplied with a volume adapter , which is divided in 2 Fix-Volumina.

By rotating the pressure knob you reach a second volume raster, which shows 50% of the total lift.

The two different volumina are named on the volume adapter and the arrow on the plunger, must be lined up with the 2<sup>nd</sup> arrow, showing on the volume adapter.

### Disassembly

**CAUTION!** The dispenser must be completely empty.

1. Screw of the volume adapter (1) of the valve head (2).
2. You now have the pressure knob (function (3)) volume adapter with the spring (4) in your hand. The PTFE-plunger is fixed to the function button (5).
3. The glass cylinder (6) is arrested into the valve head, which you can easily clean now.
4. For assembly just add all components back together again in reversed sequence

### Caution!

The upper part must be tightly screwed on the valve head, otherwise it can leak.

**Warning:** Avoid personal injury from chemicals. The instruments, suction and ejection tube may contain reagent. Always face them away from your body. Wear eye protection and protective clothing!

### Please note:

Never press the plunger downwards with force. Never raise or depress it too fast. Always leave the plunger down and empty after use.

## Cleaning and maintenance

As you can see on the sectional drawing, MINISPENSOR can be disassembled in seconds. Therefore the dispenser is optimal to clean and, if necessary, easy to change spare parts.

To ensure a proper performance and a long life-span, the dispenser should be cleaned in the following cases:

- When changing reagents
- Before storing it over a long period of time
- Daily after you have dispensed:
  - \* Substances which form crystals
  - \* Alkaline solutions
  - \* Organic solutions
  - \* anorganic oxidizable solutions ( fe. Biuret-reagent).

## Preparation for ceaning (Disassembly)

1. Place the complete dispenser (with bottle) in a sink.
2. Unscrew the apparatus from the bottle so that the suction tube will not be immersed in the liquid.
3. Tap the tube carefully against the inner side of the bottle. This allows the remaining reagent to disperse.
4. Remove dispenser from the bottle and rinse with distilled water.
5. It is now possible to disassemble the dispenser. Please refer to para „Dissasemble“

**Warning:** Avoid personal injury from chemicals. The instruments, suction and ejection tube may contain reagent. Always face them away from your body. Wear eye protection and protective clothing!

## Prevention of clogs of valves during prolonged intervals

When the dispenser is screwed onto the bottle make sure that the valves are surrounded and covered by liquid. Once unscrewed you can maintain an easy running of the valves, if you rinse it with distilled water and/or laboratory cleaner. Finally draw alcohol through the dispenser.

From time to time in order to ensure the apparatus meets its accuracy, check-ups can be done with (2x)-distilled water on a semi-micro scales. Allow temperature dependance (1 ml water 20° C = 0,998 g)!

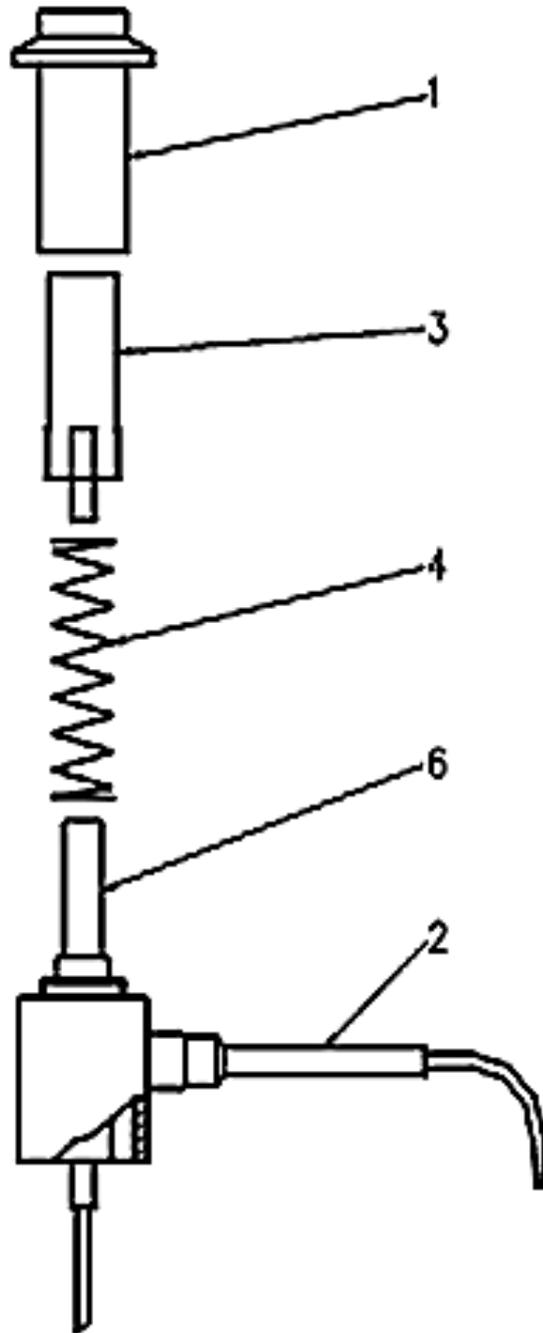
## Sterilization:

After removal of the suction tube and ejection canula the dispenser can be steam-sterilized at 121 °C and maximum 2 bar. Place the instrument in an upright position on a towel. Avoid any contact with hot metal surfaces.

## Caution:

Slowly cool down the apparatus. It can then be used again after reaching room temperature. Inspect all parts and tubing fro leakage after every sterilization. In case of deformation return the Minispensor.

It is also possible to sterilize the dispenser chemically with alcohol, formaldehyd etc.

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LABORTECHNIK GmbH