


Bedienungsanleitung / User Guide



Ansaugpumpe mit Absperrhahn / Suction pump with stopcock

Art. No. 546000

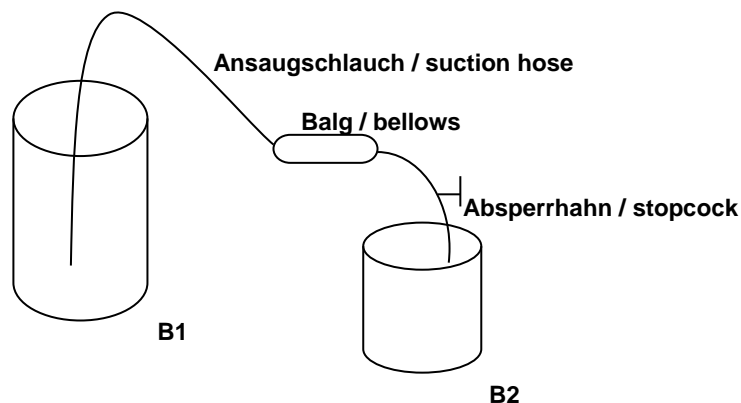

WICHTIG!

Beim Abfüllen von Flüssigkeiten muss Entnahme-Behälter B1 höher stehen als Abfüll-Behälter B2

1. Ansaugschlauch in Entnahme-Behälter halten.
2. Bei geschlossenem Hahn den Balg mit der Hand zusammendrücken und pumpen bis das Schlauchsystem komplett mit Flüssigkeit gefüllt ist.
3. Absperrhahn öffnen und Flüssigkeit in Behälter B2 fließen lassen.
4. Nach Abfüllvorgang Absperrhahn schließen.

Beim Umfüllen von Säuren, Schlauch an beiden Tüllen Seiten des Balgs mit Schlauchklemme sichern.

Suction pump with stopcock

IMPORTANT!

When filling liquids, extraction container B1 must be higher than filling container B2

1. Hold the hose into extraction container.
2. When the valve is closed hand squeeze bellows and pump up until the hose system is completely filled with fluid.
3. Open the stopcock and allow liquid to flow into tank B2.
4. After filling process close the stopcock.

When transferring acids, secure hose spouts on both side of the bellows with hose clamps.