WiseTis[®] HG-15D

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Contents

1 Introduction	8
2 Safety Instructions	8
2.1 Description of symbols.....	8
2.2 Instructions	8
3 Product description	10
3.1 Key features	10
3.2 General survey.....	10
4 Packing contents	11
5 Positioning the device and commissioning	11
6 Job description	12
6.1 Set speed.....	12
6.2 Store speed values.....	12
7 Maintenance and cleaning	12
8 Technical data	12
9 Certificate	13

1 Introduction

Thank you for purchasing our **WiseTis® HG-15D Homogenizer**.

You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**Risk of injury**

Avoid injuries of your hands by following these instructions.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Reparations and amendments should only be made by **witeg** authorized people. Contraventions exclude the warranty and may cause fatal injuries.

- Do not use this device in areas with potentially explosive atmospheres.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective clothing during the usage.
- Pull out the power cable if you don't want to use the device for a longer time.
- Dry run may destroy the dispersing tool.

- It is important that the device stands up on a horizontal and firm subsurface to avoid vibrations.
- Make sure that the device can not slip away.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superior to set the device in a good illuminated place to avoid accidents engendered by handling errors.

- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not put your hands into the device before the motor has stopped.
- Turn the device off before change dispersing tools.
- Demount dispersing tools before cleaning.

3 Product description


3.1 Key features

The **Wisdom WiseTis® Homogenizer HG-15D** is designed for nearly all media and dispersing processes in chemistry, biology, medical and pharmaceutical science.

Key features are as follows:

- Simplest operating.
- Fast commissioning.
- Long life by simply and solid construction.
- Over temperature and over current protection.
- Illuminated LCD for speed and power bar graph.
- Speed adjustable from 2000 up to 27.000 rotations per minute in steps to 50 rpm.
- The dispersing tools are made of stainless steel and PTFE and are suited for all sterilization methods.

3.2 General survey

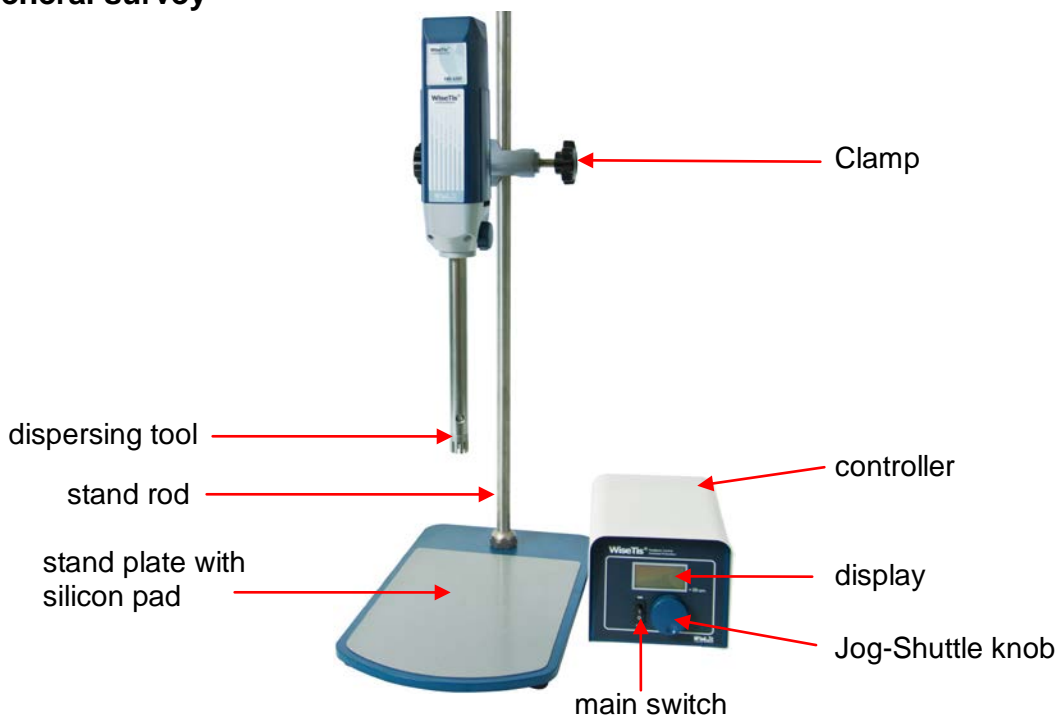4 Packing contents

quantity	denomination		
	DH.WHG02019	DH.WHG02020	DH.WHG02018
1	WiseTis® Homogenizer HG-15D (with remote controller)		
1	Allen key		
1	body holder		
1	operating manual		
1	ST110 standard plate stand		
1	CL200 clamp		
1	HT1018 dispersing tool	HT1025 dispersing tool	

5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package and perhaps soiling carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Attach the clamp to the tripod rod.
5. Mount the body holder at the clamp.
6. Stick the device at the body holder.
7. Tighten the bolts with the included Allen key carefully.
8. Stick the dispersing tool into the device and tighten the screw.
9. Connect the remote controller with both connection cables.
10. Check that each connection fits close.
11. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
12. Make sure that you follow all safety instructions.
13. Connect the device with the power outlet.
14. Switch the device on.
15. Set your values and start the process (look at job description).

Never run the device without any load and dip it into the medium which should be homogenized before powering on the device. Overspeed and resulting vibrations can damage the device.

Don't push the homogenizer constantly against the vessel, if you are homogenizing inside a glass vessel. High frequent vibrations could damage your glass vessel.

6 Job description

6.1 Set speed

- Switch the device with the main switch on.
- There was a value stored before if in the upper line turns up a value > 0. Turn the Jog-Shuttle to adjust the speed and push it to start the dispersing process.
- If there turns up "0" in the upper line the dispersing process will start by turning the Jog-Shuttle knob.
- Push the Jog-Shuttle for start and stop.

Multiply the showed speed value with 10 to get the real speed.

6.2 Store speed values

If you work with the same speed value often you can store this value and save precious time.

- Set your speed by turning the Jog-Shuttle.
- Hold the Jog-Shuttle pushed till the upper value starts blinking.
- Delete values by over-storing them with "0".
- After switching the device on you only need to push the Jog-Shuttle to start the dispersing process.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device and pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

8 Technical data

		HG-15D
Dimension (mm) (wxdxh)		60x70x215
Speed (U/min)		2000 - 27.000
Dispersing capacity		up to 2,5 liters
Motor rating input /output (W)		300 / 160
Safety device		overheat protection
Material	Housing	powder coated steel
	tools	stainless steel and PTFE (bearing)
Weight (kg)		2,5
Usable tools		HT1008, HT1010, HT1018, HT1025
Power supply		AC 230 V, 1N~, 50/60 Hz

9 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	WiseTis[®] HG-15D Homogenizer
Typ	
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de