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USER MANUAL ULTRASONIC CLEANER

PLEASE NOTE!

Drawings in this manual are for illustration purposes only and in some details may differ from the actual machine. The original operation manual is in German. Other language versions are translations from German.

INTRODUCTION

Thank you for choosing our products. To avoid injury from fire or electric shock, please ensure compliance with safety instructions when using this device. Please carefully read the instructions and make sure to have understood it well. Keep the manual near the equipment to be able to read it at any time. Always use current sources connected to the ground and providing the necessary voltage (indicated on the label on the device). If you have any doubt, let an electrician check that your outlet is properly grounded. Never use a damaged power cable. Do not open the unit in damp or wet environment, or if you are wet yourself. Protect the unit of solar radiation. Use the device in a protected location to avoid damaging the equipment or endanger others. Make sure the device is able to cool and avoid placing it too close to other devices that produce heat. Before cleaning, disconnect it. Use a soft damp cloth for cleaning. Avoid using detergent and make sure that no liquid enters the unit. No internal element of this device needs to be maintained by the user. An opening of the device without our approval leads to a loss of warranty.

GENERAL INFORMATION

Principle of ultrasonic cleaning

Under the bottom of the basin are installed ultrasonic transducers that convert electrical energy into mechanical vibrations. This leads to the formation of small bubbles which implode (cavitation). The principle of cavitation removes dirt thoroughly. It is the so called "electronic brush".

General safety tips

Only those who have read the manual and adults are allowed to use this device.

- Do not dive parts of the body or parts of living things (animals, plants) into the basin, especially when the unit is in operation.
- Avoid the contact with eyes or skin of the detergents used.
- Do not inhale the detergent.
- Do not use the device if it's empty.
- The tank of the device must be filled with minimum level of the cleaning liquid in accordance with the below table:

Modell	MINIMUM VOLUME OF THE CLEANING LIQUID
0.7M	2/3 of the tank volume
2.0M	2/3 of the tank volume
3.0M	2/3 of the tank volume
0.7WH	2/3 of the tank volume
4.0M	2/3 of the tank volume
6.0M	2/3 of the tank volume
10.0M	2/3 of the tank volume
28.0M	2/3 of the tank volume
2.0M ECO	2/3 of the tank volume
3.0M ECO	2/3 of the tank volume
6.0M ECO	2/3 of the tank volume
10.0M ECO	2/3 of the tank volume
28.0M ECO	2/3 of the tank volume
0.7	2/3 of the tank volume
2.0	1/2 of the tank volume
3.0	2/3 of the tank volume
4.0	2/3 of the tank volume
6.0	2/3 of the tank volume
10.0	1/2 of the tank volume
28.0	1/3 of the tank volume
0.7D	2/3 of the tank volume
2.0DS	2/3 of the tank volume
3.0DSP	2/3 of the tank volume
2.0ECO	2/3 of the tank volume
3.0ECO	2/3 of the tank volume

6.0ECO	2/3 of the tank volume
10.0ECO	2/3 of the tank volume
20.8ECO	2/3 of the tank volume

Other possible uses: CDs, DVDs, metal objects such as cutlery, coins, etc..

Caution

- Do not clean porous stones or pearls in the device.
- The watches must be waterproof (at least 30m).
- It is possible that loose parts of the object while cleaning.
- Objects which easily lose their color should not be cleaned in the unit.

- Do not fill the tank with water above 50 ° C.
- Do not place items that have been in contact with acids, chloride ions and other aggressive substances in the basin.
- Do not use flammable liquids or other chemicals containing chloride ions, or leading to their release (such as some disinfectants, detergents or dishwashing liquids).
- Check when unpacking the unit that it suffered no damage during the transport. If so, please contact us immediately. Keep in all cases the original packaging.
- Place the unit in a stable position on a horizontal, clean and smooth surface.
- Maintain the unit in a clean room.
- During the heating of the cleaning liquid, be sure to stir the liquid every 15 minutes so as to avoid overheating and boiling.
- During ultrasonic cleaning, you will hear the sounds of cavitation. When using the device for a long time, please wear hearing protection within a radius of 2m around the device.
- The cleaning fluid heats up when it is used long: watch the temperature if you clean objects sensitive to heat.
- Only carry the device when it is empty!

Usability

- Jewelry**
 - Necklaces
 - Rings
 - Earrings
 - Waterproof watches
- Glasses**
 - Sunglasses
 - Watchbands
- Beauty care equipment**
 - Razor blades
 - Dentures
 - Toothbrushes
 - Combs

Model	Pro-clean 0.7	Pro-clean 2.0	Pro-clean 3.0	Pro-clean 4.0	Pro-clean 6.0	Pro-clean 10.0	Pro-clean 28.0	Pro-clean 0.7D	Pro-clean 2.0DS	Pro-clean 3.0 DSP	Pro-clean 2.0 ECO	Pro-clean 3.0 ECO	Pro-clean 6.0 ECO	Pro-clean 10.0 ECO	Pro-clean 28.0 ECO
Series	memory quick							LCD Degas	Degas Sweep	Multi-function	memory quick eco				
Capacity (L)	0.7	2	3	4	6	10	28	0.7	2	3	2	3	6	10	28
Internal dimensions of tank (mm)	150x 86x 65	150x 138x 100	238x 138x 100	302x 152x 100	302x 152x 15	302x 239x 150	502x 298x 200	150x 86x 65	150x 138x 100	238x 138x 100	150x 138x 100	238x 138x 100	302x 152x 150	302x 239x 150	502x 298x 200
Power supply	230~V / 50 Hz														
Sound frequency (kHz)	40	40	40	40	40	40	40	40	37	37	40	40	40	40	40
Ultrasonic power (W)	60	60	120	120	180	240	480	60	80	160	60	80	2x80	180	6x80
Heating	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes
Drain tap	No	No	No	No	No	Yes	Yes	No	No	No	No	No	No	Yes	Yes
Handles	No	No	Yes	Yes	Yes	Yes	Yes	No	No	Yes	No	Yes	Yes	Yes	Yes
Lid	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes
„Degas“	No	No	No	No	No	No	No	Yes	Yes	Yes	No	No	No	No	No
„Sweep“	No	No	No	No	No	No	No	No	Yes	Yes	No	No	No	No	No
„Pulse“	No	No	No	No	No	No	No	No	No	Yes	No	No	No	No	No
„Power set“	No	No	No	No	No	No	No	No	No	Yes	No	No	No	No	No

Model	Pro-clean 0.7M	Pro-clean 2.0M	Pro-clean 3.0M	Pro-clean 4.0M	Pro-clean 6.0M	Pro-clean 10.0M	Pro-clean 28.0M	Pro-clean 2.0M ECO	Pro-clean 3.0M ECO	Pro-clean 6.0M ECO	Pro-clean 10.0M ECO	Pro-clean 28.0M ECO	Pro-clean 0.7WH
Series	Basic							Basic Eco					Basic Without Heater
Capacity (L)	0.7	2	3	4	6	10	28	2	3	6	10	28	0.7
Internal dimensions of tank (mm)	152x 87x 65	150x 137x 100	238x 138x 100	302x 152x 100	302x 152x 150	302x 239x 150	502x 298x 200	150x 138x 100	238x 138x 100	202x 152x 150	302x 239x 150	502x 298x 200	150x 86x 65
Power supply	230~V / 50 Hz												
Sound frequency (kHz)	40	40	40	40	40	40	40	40	40	40	40	40	40
Ultrasonic power (W)	60	60	120	2x 60	3x60	4x60	8x60	60	80	2x80	3x60	6x80	60
Heating	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	No
Drain tap	No	No	No	No	No	Yes	Yes	No	No	No	Yes	Yes	No
Handles	No	No	Yes	Yes	Yes	Yes	Yes	No	Yes	Yes	Yes	Yes	No
Lid	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes
„Degas“	No	No	No	No	No	No	No	No	No	No	No	No	No
„Sweep“	No	No	No	No	No	No	No	No	No	No	No	No	No
„Pulse“	No	No	No	No	No	No	No	No	No	No	No	No	No
„Power set“	No	No	No	No	No	No	No	No	No	No	No	No	No

GENERAL USAGE AND CLEANING INSTRUCTIONS

Cleaning tips

Simple cleaning

Just put a few drops of liquid soap in water. You can use any liquid cleaner that is not acid based.

Complete cleaning

Use proprietary cleaners to remove rust or other intense stains.

Caution

The use of liquid acid-based cleaners or alkali can lead to the formation of rust on the basin walls and even lead to the formation of holes in the walls of the unit. Be sure to use products at low pH.

PREPARATION

1. Before using the device ensure that the power supply complies with the technical requirements on the device label.
2. Place the device on a flat, even surface. Fill the tank with water ensuring that the objects to be cleaned are completely submerged. Place the objects into the basket provided.

Warning

Fill the tank up to the 2/3 mark with water. Never operate the device when empty or with insufficient water! This might damage the device.

BASIC / BASIC ECO


1. Heating ON/OFF; heating setting
2. Indicator for heating in process

3. Timer ON/OFF; Timer setting
4. Indicator for ultrasonic cleaning in process

OPERATION

1. Once the switch is rotated clockwise, the machine will enter into the time operating mode. The time can be freely modified in accordance with the indication at the edge of rotary switch on the panel.
2. The green work indication lamp will turn on and the ultrasonic will send out the working sound of "squeak" and the machine will start working.
3. The cleaning machine will turn off automatically when time is up.
4. Rotate the switch anticlockwise to start long-term operating mode. The infinity mark ∞ is indicated in the panel.
5. Green working indicator lamp is on and the ultrasonic sends out the working sound of "squeak" and starts working and will continue working.
6. When cleaning is completed, in order to power off please rotate the switch clockwise to the "off" manually.
7. Rotate the heating switch clockwise to start heating the cleaning fluid. The heating temperature can be selected in accordance with the indication at the edge of rotary switch on the panel.
8. The green heating indicator lamp will turn on and the cleaning fluid will be heated.
9. When the temperature reaches the set value, the green heating indicator lamp will turn off and heating will stop.
10. When the temperature of the fluid will be lower than set value on the rotary switch on the panel, then device will start heating again. This way, the device will maintain the set temperature.
11. To stop heating, the rotary switch must be turned to OFF position anticlockwise.

MEMORY QUICK / MEMORY QUICK ECO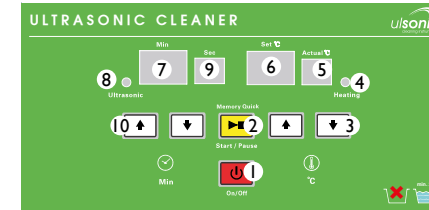

1. On/Off button
2. START/PAUSE & Memory button
3. Temperature control buttons
4. Heating process LED
5. Current temperature indicator for the cleaning liquid
6. Set temperature indicator for the cleaning liquid
7. Time indicator: displays the set time or remaining time in minutes
8. Cleaning process LED
9. Remaining cleaning time counter
10. Cleaning time adjustment buttons


USE

1. All buttons are tactile. A simple touch is enough. Press the button No. 1 to turn on. The cleaning process begins. You can now set the desired temperature and time.
2. Temperature set: Display No 5 shows the temperature of the cleaning liquid. To heat the liquid, use arrows No. 3 (up to raise the temperature, down to lower it). If the temperature is higher than the current temperature, the unit starts to heat up and the LED lights up. Once the desired temperature is reached, the diode turns off. To terminate the heating process, adjust the current temperature setting to '00'.
3. Time set: When the device is on, you see on display number 7 the time set in the previous cleaning. Press No 10 to set the time (up to increase, down to lower)
4. After setting the time, press START (No. 2) to begin the cleaning. A red LED turns on. When the countdown begins (No. 9), the cleanup is almost

complete. The cleaning will stop automatically at the end of programmed time. The time programmed by the last use appears (No. 7).

5. Pause function and memory: Press the green button Start / Pause (No. 2) to activate the memory function. When the time is up and you touch the green button, the programmed time is saved. If the cleaning process is underway and you press the green button, the cleaning is paused. To restart, repress the green button.
6. When cleaning is finished, press the ON / OFF (No. 1) button to turn off the device. Remember to unplug it after use. Empty the basin so that the unit is completely empty.

LCD DEGAS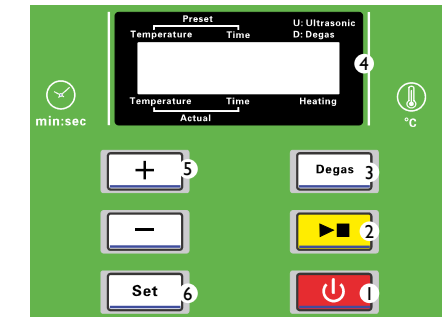

1. On/Off button
2. START/PAUSE & Memory button
3. Degas button
4. LCD display
5. Up/Down control buttons
6. Set button


OPERATION

1. All the buttons are touch buttons which only require a light touch. To activate the device and start the cleaning process, touch the power button (#1) at the front of the device. You can then select the required temperature and time settings.
2. Setting the temperature: The current temperature is shown on the LCD display. The display will flash when you press the Set button (#6) and you can now set the required temperature by pressing the

Up/Down control buttons (#5). Touch the Up or Down buttons to increase or lower the temperature setting as desired. If the set temperature is higher than the actual temperature the device will start to heat up immediately. Once the required temperature is reached, it will be maintained at this level. To terminate the heating process, adjust the current temperature setting to '00'.

- Setting the time: Press the Set button (#6). The display will start flashing. You can now program the cleaning time by pressing the Up/Down control buttons (#5) to increase or decrease the time.
- After setting the time press the Start button (#2) to initiate the cleaning process. The display shows 'U' while cleaning is in progress. Once the cleaning time has expired, cleaning will automatically stop. The display will now show the last saved time setting (Memory feature).
- Memory&Pause: Touch the green Start/Pause button (#2) to activate the Memory feature. When cleaning is complete and the green button is pressed the previous time setting is saved so that the time does not have to be re-set every time. If the green button is pressed while cleaning is still in progress, cleaning will be paused. To restart the cleaning process, press the green button again.
- Degassing: During this process ultrasound is used to expel gases (e. g. carbon dioxide) from the liquid. The number of cavitation bubbles that are required for efficient cleaning can only be formed in degassed liquids. To activate this feature press the Set button (#6). The display will show 'D'. You need to press the Set button again to stop degassing. The display will then show 'U'.
- Once cleaning is complete, switch the device off (power button #1). Remember to unplug the device after use. Empty any liquids from the tank and remove any objects.

LCD DEGAS SWEEP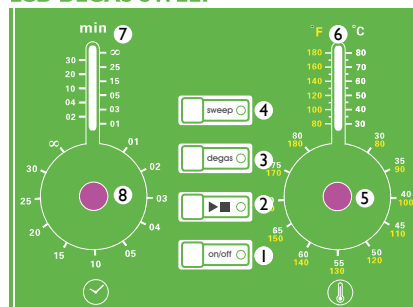

- On/Off button
- START/PAUSE & Memory button
- Degas button
- Sweep button
- Temperature control knob
- Temperature indicator
- Remaining cleaning time indicator
- Timer control knob

OPERATION


- All the buttons are touch buttons which only require a light touch. To activate the device and start the cleaning process, touch the power button (#1) at the front of the device. You can then select the required temperature and time settings.
- Setting the temperature: If the device is switched on, a flashing LED indicates the actual temperature of the cleaning liquid. To set the desired temperature, turn the temperature control knob (#5) clockwise. A green LED indicator illuminates while the device is heating up. Once the set temperature is reached it will be maintained until the device is switched off. To switch off the heating function, turn the control knob anti-clockwise to the zero position.
- Setting the time: Turn the Timer control knob (#8) to set the desired cleaning time. Pressing the Start button (#2) will activate the device and the cleaning process will start. A flashing LED indicates the remaining cleaning time. At the end of the cleaning time the cleaning process will stop automatically. The display will now show the last saved time setting (Memory feature).

- Pause: If you press the Pause button (#2) the cleaning process will be interrupted until the Pause button is pressed again.
- Degassing: During this process ultrasound is used to expel gases (e. g. carbon dioxide) from the liquid. The number of cavitation bubbles that are required for efficient cleaning can only be formed in degassed liquids. To activate this feature press the Degas button (#3). Press the button again to stop degassing.
- Sweep: This optional Sweep feature optimises the sound field distribution in the cleaning liquid through frequency modulation. Pressing the Sweep button (#4) will activate the feature. To turn the feature off, press the button again.
- Once cleaning is complete, switch the device off (power button #1). Remember to unplug the device after use. Empty any liquids from the tank and remove any objects.

OPERATION

- All the buttons are touch buttons which only require a light touch. To activate the device and start the cleaning process, touch the power button (#1) at the front of the device. You can then select the required temperature and time settings.
- Setting the temperature: If the device is switched on, the actual temperature of the cleaning liquid is indicated on the display. Turn the control knob (#14) to set the required temperature. The heating process will start automatically and the display will indicate that heating is in progress. The set temperature is maintained as long as the device is switched on. To stop the heating, turn the Temperature control knob anti-clockwise to the zero position.
- Setting the time: Turn the Timer control knob (#7) to set the desired cleaning time. Pressing the Start button (#2) will activate the device and the cleaning process will start. The remaining time is shown on the display. At the end of the cleaning time the cleaning process will stop automatically. The time display now shows the last saved time setting (Memory feature).
- Pause: If you press the Pause button (#2) the cleaning process will be interrupted until the Pause button is pressed again.
- Degassing: During this process ultrasound is used to expel gases (e. g. carbon dioxide) from the liquid. The number of cavitation bubbles that are required for efficient cleaning can only be formed in degassed liquids. To activate this feature press the Degas button (#5). The corresponding LED will illuminate. To disable degassing press the button again. The corresponding LED will go off.
- Sweep: This optional Sweep feature optimises the sound field distribution in the cleaning liquid through modulation of frequencies. Press the Sweep button (#4) to enable this feature. The corresponding LED will light up. To disable the feature press the Sweep button again. The corresponding LED will go off.

MULTIFUNCTION

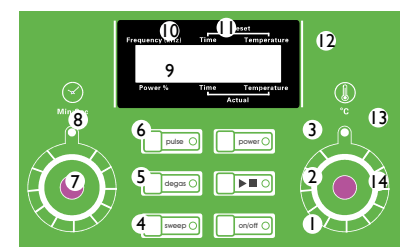- On/Off button
- START/PAUSE & Memory button
- Ultrasonic Power setting button
- Sweep button
- Degas button
- Pulse button (wave frequency setting)
- Timer control knob (cleaning time setting)
- Cleaning process LED
- Power indicator
- Cleaning frequency indicator
- Cleaning time indicator
- Temperature indicator for the cleaning liquid
- Heating process LED
- Temperature control knob (cleaning liquid temperature setting)

7. Setting the wave frequency: Pressing the Pulse button (#6) increases the wave frequency to 120% of the default value. The corresponding LED light will illuminate. Pressing the Pulse button again will reduce the value to 100%. The corresponding LED will go off.
8. Setting the ultrasonic power: The default value is set to 40% and can be increased to 70% or 100%. To change the value, press the Power button (#9) until the desired value is displayed.
9. Once cleaning is complete, switch the device off (power button #1). Remember to unplug the device after use. Empty any liquids from the tank and remove any objects.

BASIC WITHOUT HEATER

OPERATION

1. Once the switch is rotated clockwise, the machine will enter into the time operating mode. The time can be freely modified in accordance with the indication at the edge of rotary switch on the panel.
2. The green work indication lamp will turn on and the ultrasonic will send out the working sound of “squeak” and the machine will start working.
3. The cleaning machine will turn off automatically when time is up.
4. Rotate the switch anticlockwise to start long-term operating mode. The infinity mark ∞ is indicated in the panel.
5. Green working indicator lamp is on and the ultrasonic sends out the working sound of “squeak” and starts working and will continue working.
6. When cleaning is completed, in order to power off please rotate the switch clockwise to the “off” manually.

REGULAR CONTROL OF THE DEVICE

Check regularly the device to see if it presents any damages. If there is any damage, please stop using the device. Please contact your customer service to solve the problem.

What to do in case of a problem?

Please contact your customer service and prepare the following information:

- Invoice number and serial number (the latter is to be found on the technical plate on the device).
- If relevant, take a picture of the damaged, broken or defective part.
- It will be easier for your customer service clerk to determine the source of the problem if you give a detailed and precise description of the matter. Avoid ambiguous formulations like “the device does not heat”, which may mean either that it does not heat enough or that it does not heat at all. The more detailed your information, the better the customer service will be able to solve your problem rapidly and efficiently!

CAUTION: Never open the device without the authorization of your customer service. This can lead to a loss of warranty!

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