ISO 9001
CERTIFIED

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Instruction Manual

Please read before use!

pipetus®-standard

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electrical pipetting aid

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Contents

	<i>Page</i>
<i>1. Safety Instruction</i>	<i>3</i>
<i>2. Operating principle</i>	<i>6</i>
<i>3. Use</i>	<i>7</i>
<i>4. Pipetting</i>	<i>8</i>
<i>5. Maintenance</i>	<i>9</i>
<i>6. Available spare parts</i>	<i>10</i>

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1. Safety Instructions

Attention: Before use please read the safety instructions and observe them carefully. Before use every member of staff must be familiar with these safety guidelines and they must be kept available at all times. This instruction manual cannot, however, cover every possible safety hazard. It is every user's responsibility to observe all health and safety precautions and to determine and observe restrictions before use of the unit.

- Observe all general safety practices and guidelines for laboratories, such as the use of protective clothing, safety glasses and gloves when handling hazardous chemicals.

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Attention: If the unit is opened, the warranty is invalidated and any resultant claims void.

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- *The filter supplied contains a hydrophobic 0,2 µm membrane for air filtration in either direction.*
- *Only have the unit professionally serviced.*
- *Never use undue force!*

Attention: *Only ever use the original accessories and replacement parts to avoid damaging the unit.*

2. Operation principle

pipetus®-standard is an electrical pipetting aid for all conventional pipettes from 0,1 up to 100 ml. pipetus®-standard is controlled by separate buttons to fill (IN) and dispense

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(OUT). The speed of filling or dispensing can be regulated by pressure to the respective button. For working with pipettes EX (TD = to deliver) disconnect the green hose from the pump to avoid that the pipette is blown out by the pumps pressure. Like that the pipette runs out by gravity.

3. Use

Connect pump to a 220 voltage resp. 110 outlet.

Caution: Check voltage between pump and power supply.

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4. Pipetting

Insert the pipette into the pipetholder to ensure a tight fit. When using EX (TD = to deliver) adjusted pipettes, the green rubber tube has to be disconnected from the pump. When using blow out pipettes, make sure that the green tube is fixed to the corresponding outlet at the pump. For filling press the IN button, for emptying press the OUT button. Finger pressure on the button controls the working speed. Additionally the pump's power can be regulated with a nut (part no. 990 42 0 1).

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5. Maintenance

When normally using the pipetus®-standard, it is maintenance free. Depending on the use, the filter will require replacement. The nose piece assembly (990 37 00), except the filter (990 37 04), can be sterilized by autoclaving at 121°C.

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6. Available spare parts

990 37 00	pipetholder complete
990 37 01	pipetholder case
990 37 04	round filter set (with one-sided connection, 5 filters)
990 32 01	silicon adapter
990 32 02	safety valve assembly
990 31 00	handle complete
990 31 06	hose connector
990 41 04	microfibre-filter set for pump (5 filters)
990 42 01	nut M 5
990 42 03	rubber foot
990 34 00	pump without filter 230 V
990 44 00	pump with filter 230 V
990 54 00	pump without filter 120 V
990 64 00	pump with filter 120 V