made in germany  made by ritter

# Ripette®

DISPENSER

liquid handling

Artikel-Nr. 42001-0000

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Operating Manual

Mode d'emploi

Manual de instrucciones

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# Ripette®

You have made a good choice with the Ripette®. This dispenser fits optimally in your hand and is ideally suited for serial dosing due to its light weight.

With the classic 8 tip sizes, you can flexibly select by 5 stroke adjustments your desired dosing volume from 40 dosing programs. You can view all settings and volumes at a single glance through the table on the Ripette®.


Although the Ripette® is intended to be ideally used with Ritips® and Ritips® Professional, you can still work accurately and compatibly with PD-Tips® from Brand, Combitips® and Combitips® Plus from Eppendorf™.

The Ripette® is made from high-quality materials and is highly resistant. Thanks to a purely mechanical, maintenance-low and robust construction, the Ripette® has proven itself on the market for years as a durable product. Nevertheless, we recommend a calibration every 12 months in order to assure the correctness of the dispensed volumes.

The Ripette® is a dosing device for operator-controlled or defined dispensing of liquids. Using Ritips® and Ritips® Professional (referred to below as "tips"), volumes from 1 µl to 5000 µl can be dispensed repeatedly with a high degree of accuracy.

## Contents

Operation .....	3
I. Dosing Table and Volume Selection	
II. Operating Elements	
III. Inserting the Tip	
IV. Filling the Tip	
V. Dosing	
VI. Removing the Tip	
Troubleshooting .....	6
Maintenance and Care .....	7
Guarantee .....	7
Safety Regulations .....	8
Technical Data .....	9
Ordering Information .....	10

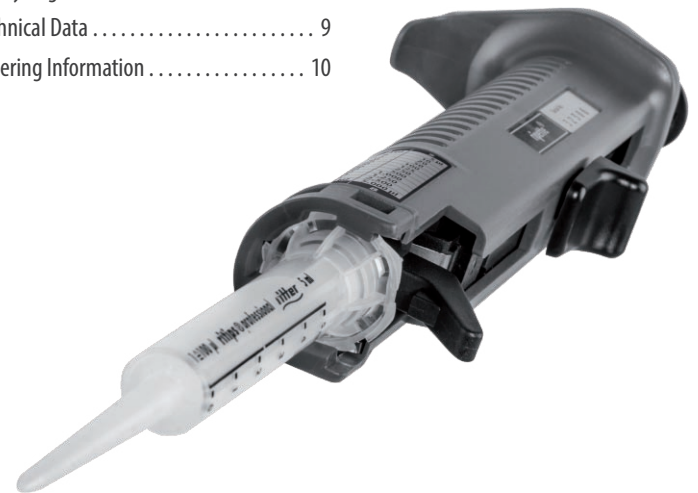# Operation

NO	1	2	3	4	5
0.05 ml	1 µl	2 µl	3 µl	4 µl	5 µl
0.10 ml	2 µl	4 µl	6 µl	8 µl	10 µl
0.20 ml	4 µl	8 µl	12 µl	16 µl	20 µl
0.50 ml	10 µl	20 µl	30 µl	40 µl	50 µl
1.00 ml	20 µl	40 µl	60 µl	80 µl	100 µl
1.25 ml	25 µl	50 µl	75 µl	100 µl	125 µl
2.50 ml	50 µl	100 µl	150 µl	200 µl	250 µl
5.00 ml	100 µl	200 µl	300 µl	400 µl	500 µl
10.00 ml	200 µl	400 µl	600 µl	800 µl	1000 µl
12.50 ml	250 µl	500 µl	750 µl	1000 µl	1250 µl
25.00 ml	500 µl	1000 µl	1500 µl	2000 µl	2500 µl
50.00 ml	1000 µl	2000 µl	3000 µl	4000 µl	5000 µl
STEPS	48	23	15	11	8

## I. Dosing Table and Volume Selection


1. First of all, use the table (3) on the Ripette® to look for the requested dosing volume in µl.
2. The topmost line (NO) corresponds to the positions 1 to 5 of the adjusting wheel (2).
3. For the suitable tip size (7), please refer to the left-hand column.
4. For the maximum quantity of dispensing steps, please refer to the bottom most line (STEPS).

Note:

- The dosing table of the manual dispenser conveniently displays the total dispensing volume proportional to the dispensing step and tip sizes.

Example as represented:

- The dispenser dispenses in program number 4 (NO) exactly 200 µl in 11 dosing steps (STEPS) when a 2.5 ml tip is used. NO 1 corresponds to 50 µl in maximum 48 steps with the same tip size or 500 µl when a 25 ml tip is used.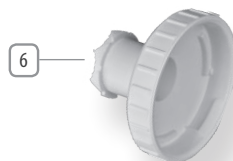

## II. Operating Elements

**Dosing button (1)** By pressing the dosing button down to the stop, the adjusted volume is dispensed.

**Adjusting wheel (2)** By moving the adjusting wheel into the position 1 to 5, the requested dosing volume is set in compliance with table 1.

**Dosing table (3)**

**Filling slide (4)** By pushing the slide upwards, the inserted tip is filled up.



**Stopping lever (5)** The stopping lever is used to secure the inserted tip.

## III. Inserting the Tip

1. Push the filling slide (4) right down to the stop.
2. Push the stopping lever (5) upwards.
3. Exert a slight pressure to insert the tip (7) which has to lock into place.
4. Push the filling slide down a bit and press down the stopping lever in position (d) so that it clicks into place.
5. Use the adjusting wheel (2) to set the required dosing volume.

Note:

- Dispenser tips 25 and 50 ml can only be inserted/locked into place by means of the enclosed adapter in the Ripette®.
- Place the adapter on the tip and fasten by turning (bayonet socket). After use of the tip, the adapter can be removed again and be reused as often as desired and even be sterilized, if required.
- A 0.05 ml tip may be used only with a 10 µl pipetting tip in its place.

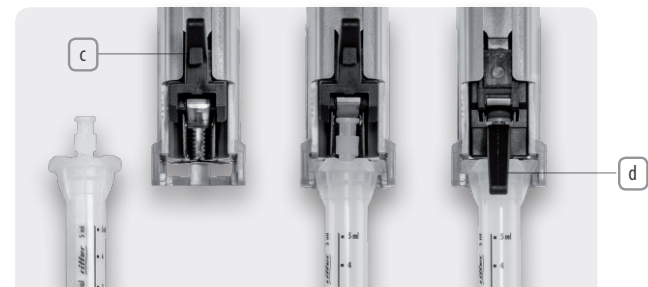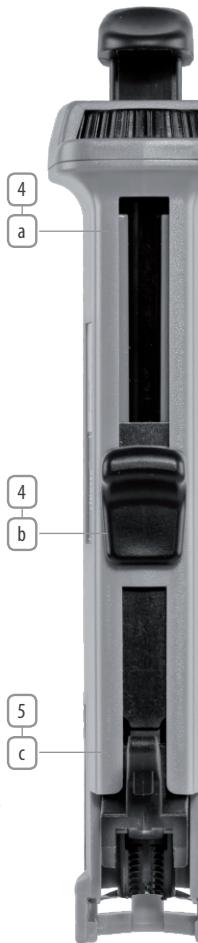## Operation

### IV. Filling the tip

1. Immerse the tip into the liquid.
2. The tip has to be filled and emptied (moistened) by pressing the filling slide (4) into the position (a) and back to position (b), in order to achieve a humidity balance in the dead air volume.
3. Fill the tip slowly and without air bubbles.

Note:

- In order to prevent a negative pressure and a saturation of the liquid with minute air bubbles when filling the tip, the filling slide has to be pushed up slowly.
- The tip does not need to be filled up completely.
- An air bubble which might be in the upper part of the cylinder will have no influence on the accuracy.
- Tips 0.5 ml, 1.25 ml and 2.5 ml can be used with fixed yellow or transparent pipetting tips. Filling from vessels with narrow necks is thus facilitated, and the dosing of smaller volumes will be more precise.
- Due to physical reasons, the contents of the first dosing step and the last dispensing step (residual stroke lock) must be discarded.

### V. Dosing

1. Check the volume setting at the adjusting wheel (2).
2. Wipe the tip with a lint-free cellulose wadding.
3. Discard the first dispensing step.
4. Place the tip at an angle to the wall of the vessel.
5. Press the dosing button (1) down to the stop.

Note:

- The tip must be changed when the reagent is changed.
- Discard the residual liquid by pressing the filling slide (4) down slowly to the stop (b).

### VI. Removing the Tip

1. Empty the tip completely.
2. Push the stopping lever (5) upwards and remove the tip carefully.

## Troubleshooting

Problem	Cause	Solution
The tip cannot be inserted.	The filling lever (4) is not pushed all the way down.	Push the filling lever into position (b) until it stops.
The stop lever (5) cannot be pushed down.	The tip is not properly inserted.	Remove the tip. Push the piston of the tip all the way in. Push the filling lever into position (b).
The filling lever cannot be pushed up.	The stop lever has not clicked into place in the tip.	Press the stop lever until it stops.
The dosing volume is indeterminate.	The adjustment wheel (2) has not clicked into place.	Click adjustment wheel into place.
An air bubble in the tip is becoming larger.	The tip is leaky.	Replace the tip.
The tips 25 ml and 50 ml cannot be inserted.	The adapter (6) is not mounted on the tip.	Remove enclosed adapter from the Ripette® original packaging and attach it on the tip.

## Maintenance and Care

- The Ripette® is factory-calibrated and works maintenance-low.
- Potential clogging is removed best of all with warm water or isopropyl alcohol.
- To disinfect the Ripette®, conventional disinfectants suitable for polycarbonate can be used.
- The device should not be autoclaved.
- The device should not be disassembled.
- No technical modifications should be made.
- A calibration every 12 months is recommended to check the correctness of the dispensed volumes.

## Guarantee

- Ritter Medical issues a guarantee during 1 year from the date of purchase for all Ripette® devices with serial numbers, subject to proper handling, maintenance and care according to the instructions for use.
- If problems cannot be corrected according to the troubleshooting information indicated, please send the Ripette® to your specialist dealer for replacement or repair.

## Safety regulations

- Read the instructions for use before using the Ripette®.
- Check the condition of the Ripette® before use. In the event of any problems, follow the instructions in the "Troubleshooting" section first.
- The Ripette® is used exclusively for dosing liquids into suitable containers.
- As a basic principle, work should be performed such that the device is held pointing down, in order to not endanger neither the user nor other persons.
- Avoid contact with the opening of the tip while working with aggressive media – reagent manufacturer's information should be observed.
- General warnings and safety instructions (e.g. wearing protective clothing) should be followed. When working with infectious or hazardous samples, Good Laboratory Practice (GLP) must be maintained.
- When working with viscous solutions, the leak tightness of the tip can no longer be guaranteed following repeated use.

## Technical Data

### Weight of the Ripette®

Weight (without tip): approx. 105 g.

### Accuracy and Precision

Liquid: double distilled water.

Reference temperature: 20 °C to 25 °C, ± 0.5 °C constant.

Measurement: in accordance with German standard EN ISO 8655.

	NO	1	2	3	4	5	S. E.*	R. E.**
	STEPS	48	23	15	11	8		
Tip Size	0.05 ml	1 µl	2 µl	3 µl	4 µl	5 µl	± 1.0	± 1.5
	0.10 ml	2 µl	4 µl	6 µl	8 µl	10 µl	± 1.0	± 1.0
	0.20 ml	4 µl	8 µl	12 µl	16 µl	20 µl	± 1.0	± 1.0
	0.50 ml	10 µl	20 µl	30 µl	40 µl	50 µl	± 1.0	± 0.6
	1.00 ml	20 µl	40 µl	60 µl	80 µl	100 µl	± 0.9	± 0.4
	1.25 ml	25 µl	50 µl	75 µl	100 µl	125 µl	± 0.8	± 0.4
	2.50 ml	50 µl	100 µl	150 µl	200 µl	250 µl	± 0.7	± 0.4
	5.00 ml	100 µl	200 µl	300 µl	400 µl	500 µl	± 0.6	± 0.3
	10.00 ml	200 µl	400 µl	600 µl	800 µl	1000 µl	± 0.5	± 0.3
	12.50 ml	250 µl	500 µl	750 µl	1000 µl	1250 µl	± 0.5	± 0.3
	25.00 ml	500 µl	1000 µl	1500 µl	2000 µl	2500 µl	± 0.5	± 0.3
	50.00 ml	1000 µl	2000 µl	3000 µl	4000 µl	5000 µl	± 0.5	± 0.25
	Dosage Volumes							

\* Systematic error at 10 % of the nominal volume in %.

\*\* Random error at 10 % of the nominal volume in %.

The data for 0.05 ml tips are based on a measurement carried out with a 10 µl pipetting tip.

Tips and dispenser have received the certificate of conformity with the following symbol, in accordance with the German Weights and Measures Regulations as amended on 12<sup>th</sup> August 1988:

For questions concerning Ripette® Ritips® and Ritips® Professional, please contact:

Ritter GmbH Medical  
 Phone: +49 8232 5003-45  
 Fax: +49 8232 5003-48  
 contact@ritter-medical.de

## Ordering Information

The Ripette® is intended for use with Ritips® and Ritips® Professional. When using compatible dispenser tips from other manufacturers, the user must check suitability before use and perform function and volume testing.

The use of Ritter's dispenser tips is therefore recommended. Optimal dispenser tips can be obtained from your specialist dealer under the following item numbers:

Item Description	Sales Unit	Article no. standard	Article no. bioclean®
Ritips® 0.05 ml	100	40002-0000	40002-0019
Ritips® Professional 0.1 ml	100	40007-0000	40007-0011
Ritips® Professional 0.2 ml	100	40007-0001	40007-0012
Ritips® Professional 0.5 ml	100	40007-0003	40007-0013
Ritips® Professional 1.0 ml	100	40007-0004	40007-0014
Ritips® 1.25 ml	100	40002-0102	40002-0111
Ritips® Professional 2.5 ml	100	40007-0005	40007-0015
Ritips® Professional 5.0 ml	100	40007-0006	40007-0016
Ritips® Professional 10.0 ml	100	40007-0007	40007-0017
Ritips® 12.5 ml	100	40002-0005	40002-0014
Ritips® Professional 25.0 ml	25	40007-0009	40007-0019
Ritips® Professional 50.0 ml	25	40007-0010	40007-0020
Ripette® with adapter	1	42001-0000	
Adapter for tip sizes 25 ml and 50 ml	50	43001-0047	autoclavable

Subject to technical modifications.

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## Products and Development

Ritter GmbH was founded in 1965. The medium-sized independent family-owned company is based in Schwabmünchen, Bavaria.

At its 25,000 m<sup>2</sup> production facility, the company's approximated 280 employees develop and produce high quality plastic products for the world market, certified under ISO 9001:2008, ISO 50001:2011 and EN ISO 13485:2003 + AC:2009.

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## Ecological

Ritter GmbH only applies certified, pollution-free plastics. When applicable, biodegradable as well as recyclable plastics are used for the production of high-quality products from Ritter Medical.

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Die Ritter GmbH wurde 1965 gegründet. Das mittelständische Familienunternehmen ist in Schwabmünchen (Bayern) ansässig.
Ritter GmbH a été fondée en 1965. Cette entreprise familiale de taille moyenne est située à Schwabmünchen, en Bavière.
Ritter GmbH fue fundada en 1965. Esta empresa familiar de tamaño mediana tiene su sede en Schwabmünchen, Baviera.