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## Operating Manual

for

# TITREX 2000

## Digital Burette

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**Please study carefully and  
follow Step by step!**

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# Index

<b>1</b>	<b>Safety Instructions</b>	<b>1</b>
<b>2</b>	<b>Use</b>	<b>2</b>
<b>3</b>	<b>Commissioning of the instrument</b>	<b>3</b>
3	Installation of Discharge tube	3
3	Installation of Suction tube	3
3	Air purging	3
4	Switch on	4
<b>4</b>	<b>Using the burette</b>	<b>4</b>
4	Titration	4
5	Dosing	5
5	Draining	5
5	Charging	5
<b>6</b>	<b>Titrexsoft</b>	<b>6</b>
6	Installation manual	6
6	Preparation	6
7	Storage of data	7
7	Data transfer	7
<b>8</b>	<b>Calibration Mode</b>	<b>8</b>
8	Calibration without software and PC	8
9	Calibration with software and PC	9
<b>9</b>	<b>Common instructions to user</b>	<b>9</b>
<b>10</b>	<b>Disassembly / Assembly</b>	<b>10</b>
<b>11</b>	<b>Cleaning</b>	<b>11</b>
<b>12</b>	<b>Accessoires</b>	<b>12</b>

*Technical modifications reserved*

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## Safety Instructions

1. Every user must study these instructions carefully before using the instrument.
2. General operation and safety instructions must be followed. Eye protection, protective clothing and glove must be worn.
3. The instrument must only be used for intended purpose and utilized within the component tolerances of material resistance.
4. Always work with the instrument so as not to endanger the operator or other persons.
5. Do not apply force to the instrument.
6. Do not drop the instrument.
7. Do not immerse the instrument in liquids.
8. Only use original spares and accessories.
9. Repairs must only be carried out by trained personnel. Do not open the instrument. Such actions will nullify the guarantee!
10. Do not use the instrument in explosive environments.
11. Do not expose the instrument to temperatures exceeding 50°C. Do not autoclave !
12. Keep the instrument away from fire or flames.

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### ATTENTION

For a problemless functionality please observe a firm seat of cup nut. If necessary please redraw.

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The TITREX 2000 Digital Burette is intended to be used for titration and dosing within following limits:

15 – 40 °C of reagent and instrument

When using according to the instructions, only these parts of the Digital Burette will come into direct contact with following chemically resistant material:

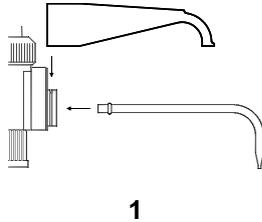
- FEP (Suction hose, Discharge tube)
- Tefzel® (Valve block)
- Borosilicate glass (Valves, KPG Cylinder)
- Ceramic, Platinum-Iridium (Valves)

Never use the instrument with

- Hydrofluoric acid
- liquids which corrode PTFE, FEP, Tefzel®, Platinum-Iridium or where catalytic changes to these material can occur
- crystallizing, particulating or precipitating solutions (e.g. suspensions)

The including adapters are made of polypropylene and shall only be used with non PP-aggressive media!

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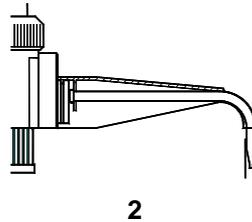
## Commisioning of the instrument Installation of Discharge Tube


Put supplied Discharge Tube into the opening and clip-on the protective sleeve (1). Fix the Discharge Tube by pressing slightly into the protective sleeve (2). Pay attention that the bulge is placed between valve body and fixing bridge of the protective sleeve (3).

## Installation of Suction tube

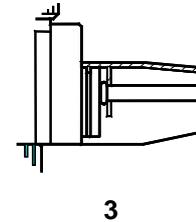Cut suction tube to desired length and put it into the smaller socket of the bottom part of the instrument. Check for firm seating.

## Air purging

Seat instrument onto the flask, and turn the plunger to below stop-point. Make short titration and dispensing moves till the medium comes out without bubbles.

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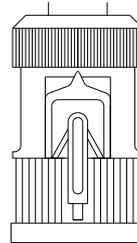
## Using the burette

### Titration

The titration mode measures the emitted volume in 0,1 ml steps.

Turn the Discharge Tube in discharge position.

Now you can titrate 99,99 ml without renewed knob-pushing.

Did you reach the 99,99 ml point, you have to set the display “00,00” before you can measure titration beyond. Push button “T” and shortly after button “Res”. Now you can go on with titration.

You can aspirate new liquid during titration as oft as necessary. If you want to aspirate certain plenty of liquid you can measure by pressing button „D“ (More information at “Dose”)

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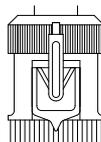
## Charging

If necessary, charge the built-in NiCd batteries with the supplied charging unit. First put the charging unit into main socket and then the plug into the bush at the back of the burette. Please let change the batteries after 3 years.

## Dosing

The dosing mode measures the aspirated volume in 0,1 ml Steps.

During dosing mode a dot is blinking between 1<sup>st</sup> and 2<sup>nd</sup> digit.

Firstly, turn the plunger completely down and set display "00,00". Push therefore button "D" and shortly after button "Res".

Now aspirate your desired volume and then turn the plunger completely down to eject the measured volume.

## Draining

To drain the discharge tube without losses turn the selector switch to position *draining*. The remaining liquid in the discharge tube flows back into the flask.

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### Features

1. When you start the first time, please insert the password "a" under admin(instrator)
2. Please select your environment data
3. You are now on the main working desk
4. Please press "service" and select your connecting parameters under "settings":
  - serial number from backside of TITREX
  - serial interface like connected: 1 – 8
  - balance: 1 (standard)
  - standard comp-value TITREX 2000 (1995 to 2003): 4130
  - Standard comp-value TITREX 2004 (after 2004): 8236
5. Please select "service", "TITREX", "comp-value" and switch on the TITREX.

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6. Please read out the comp-value to check.
7. Please click „help“ if comp-value is not readable or contact your customer service.

## Storage of data

The TITREX 2000 possess the facility to store up to 1000 values in the chip. Transfer this values with interface cord and „Titrexsoft“ to conventional PC.

Store values by pushing button „MEM“. Also stored is information if value is from titration or dosing.

Delete the actual displayed value by pushing first „MEM“ and in addition „D“ button.

Delete all values by pushing „RES“ button instead of „D“ button.

## Data transfer

You can transfer your stored values with data wire and Titrexsoft to your PC and import into a Excel table.

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Click „Save“ and select folder in which you want to store data.

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## Calibration Mode

The instrument is factory calibrated to 20°C aqueous solutions. For solutions of varying density and viscosity, the TITREX 2000 can be adjusted by the user in case of changed conditions provided a laboratory balance with a resolution of  $<1$  mg is available.

### Calibration without software and PC:

The instrument is factory calibrated to 20°C aqueous solutions. For solutions of varying density and viscosity the TITRAMAX can be adjusted by the user, provided that laboratories balance with a resolution of  $<1$  mg is available.

- Weight out 5 titrations of nominal volume (50.00 ml) and calculate the volume with temperature correction factor (order our justification procedure for titrex)
- Average the 5 resultant values
- Depress and hold button „RES“ until in the display appears „CAL“, release button and in the display the dots flashes between the first and second and as well as between the third and fourth digit
- Adjust with „D“ and „T“ the calculated average
- To activate the calibration value press button „MEM“
- Press button „RES“ to return to operation mode
- During the work with this calibration value, in the display flashes the dot between 2<sup>nd</sup> and 3<sup>rd</sup> digit
- In order to return to the factory settings, depress button „RES“ until in the display „CAL“ appears, then deactivate the calibration value by pressing the button „MEM“ and return to operation mode via pressing „RES“

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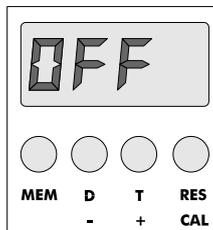
## Calibration with software and PC:

- Connect device with PC and start “Titrexsoft”. Follow “Titrexsoft” installation manual.
- Press “Service” and then “Comp-value” button.
- Switch the device on.
- Read out the comp-value.
- If this is not possible, look at „Help“ for a solution
- Calculate average mass of the 10 masses of weighed nominal volumes.
- Type in average mass in field “Balance”.
- Click “write” button in field “Calculate Comp-Value”

## Common instructions to user

In case the burette remained for longer period screwed to the flask and not in regular use, we recommend to purge the instrument prior to again titration or dosing.

Caution !! The instrument is not equipped with an „off switch“!

The TITREX 2000 automatically is switching off when you do not use it for 3 minutes. The display of the instrument indicates „OFF“ before switching out. Simply press button „D“ or button „T“ to reactivate it. In the display now appears the last titration or dosing value.

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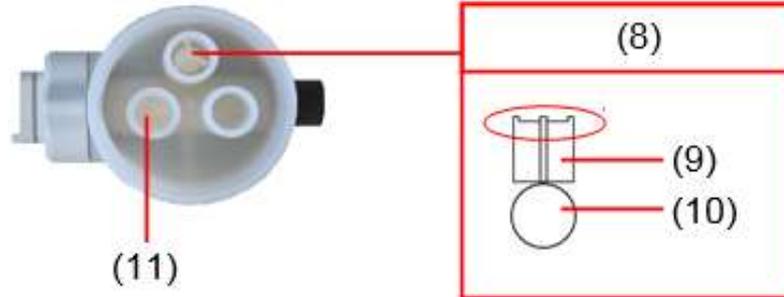
(10) Valve ball

(11) Ejection valve

Place the valve ball (10) and the valve star (9) into the suction valve (8). Check that the spikes of the valve star (9) point upwards.

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## Cleaning

The instrument should be cleaned regularly but always on

- reagent change
- long storage
- use of crystallizing media

## Warning !

**When cleaning observe all safety instructions. Components of the instrument are contaminated with titrating liquid! Suitable protective clothing, safety glasses and gloves must be worn!**

- Drain instrument and place it along with the bottle inside a washbasin
- Unscrew instrument from the bottle and pull out suction tube out of liquid
- Demount tube, let it drain and mount it again or tap carefully at the inner flask for draining
- Turn hand wheels like at air purging short up and down until there comes only a modicum of liquid
- Finally rinse TITREX with distilled water or ethanol

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witeg Labortechnik GmbH