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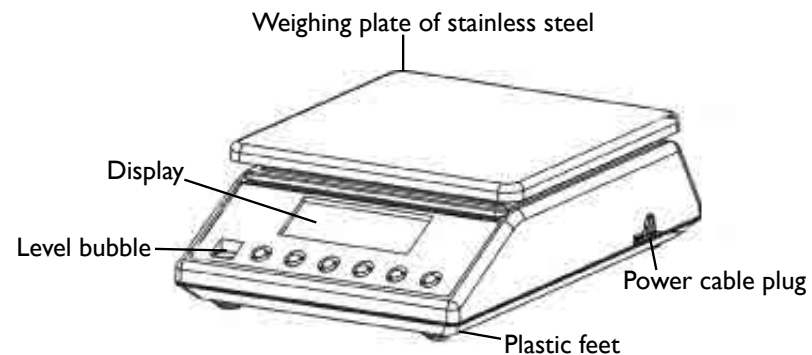
II. FUNCTIONS OF THE KEYS

1. ON/OFF

Press the ON / OFF button to start the unit. Hold the button pressed for 3 seconds to turn off the scale.

2. UNIT

The unit change is possible by pressing the MODE button. Then press up to reach the desired unit.

There are 13 units (g / ct / lb / oz / dr / gn / ozt / dwt / MM / tl.J / tl.T / tl.H / t).

3. COUNTING (PCS)

Counting mode: Press PCS to enter the counting function. The scale switches from the weighing mode to the counting mode. By pressing this button, you can switch from one mode to another.

4. TARE

This button is used to calibrate the scale. Place an empty container on the weighing plate until the balance indicates a stable value. Then press „ZERO“ Your balance now displays „0“. To exit the tare function, remove the container and repress „ZERO“: the scale is back to normal.

5. GREEN BACKLIGHT OF THE DISPLAY

Backlight of the display is turning off automatically if there is no load on the scale. The scale goes into standby mode. During weighing, the backlight of the display is turned on.

6. ENERGY SAVING FUNCTION

The scale turns off automatically after the time scheduled (2, 5 or 8 minutes of inactivity). This function is turned off by factory default settings. In order to turn it on, full calibration procedure must be performed in accordance with the below instruction. Calibration requires standard weights 5 kg and 7.5 kg / or 10 kg depending on the scale model:

1. Turn OFF the scale.
2. Press TARE and ON/OFF simultaneously for about 5 seconds. Display will show SET and then oFF (FOO).
3. Press ON/OFF – display will show 0(F01).
4. Press ON/OFF – display will show 1.0(F02).
5. Press ON/OFF – display will show 0.5(F03).
6. Press ON/OFF – display will show 1.0(F04).
7. Press ON/OFF – display will show X(F05) – where X can be 0 (AUTO-OFF is off) or 2 (scale will turn off after 2 minutes of inactivity) or 5 (scale will turn off after 5 minutes of inactivity) or 8 (scale will turn off after 8 minutes of inactivity). X can be changed by pressing MODE button.

8. Press ON/OFF – display will show 1(F06).
9. Press ON/OFF – display will show 4(F07).
10. Press ON/OFF – display will show 75000(CAP) or 100000(CAP) depending on the scale model.
11. Press ON/OFF – display will show 1(div).
12. Press ON/OFF – display will show 0.1(DEC).
13. Press ON/OFF – display will show on(LCP).
14. Press ON/OFF – display will show 5000.0(WE1).
15. Press ON/OFF – display will show 7500.0(WE2) or 10000.0 (WE2) depending on the scale model.
16. Press ON/OFF – display will show 30.....(SCL).
17. Press ON/OFF – display will show LoAd and blinking 50000 (WE1).
18. Put on the scale standard weight 5 kg – display will show some numbers and then PASS and the weight value.
19. Then the display will show LoAd and blinking 75000 (WE2) or LoAd and blinking 100000 (WE2) depending on the scale model.
20. Put on the scale standard weight 7,5 kg or 10 kg depending on the scale model – display will show some numbers and then PASS and the weight value.
21. End of calibration.

III. USE

Press the ON / OFF to turn on the scale. After turning on the scale, display first shows “000000” to „999999”, and then “0.0g”. Let the scale to acclimate for 15min to achieve optimum precision results.

OVERLOAD

When the weight of the object exceeds the capacity of the scale, it displays “HHHH”. Remove immediately the weight of the scale to prevent damages.

TARE

Place an empty container on the weighing plate until the balance indicates a stable value. Then press „ZERO“ Your balance now displays „0“. To exit the tare function, remove the container and repress „ZERO“. The scale is back to normal.

COUNTING FUNCTION

Put a number of countable parts on the scale (10, 20, 30, 40, 50, 60, 70, 80, 90, 100) and then press PCS. You pass in counting mode. Press the PCS button to select the number corresponding to the number of pieces set. Press the ON / OFF button to confirm. Your balance is now in counting mode. To return to weighing mode, simply press the MODE button.

IV. CALIBRATION (requires standard weight 5kg)

1. Hold the MODE button pressed for 4 seconds to switch to the calibration mode, the total capacity of the scale is displayed.
2. Place the corresponding standard weight on the scale until the word „PASS“ is displayed.

Your scale is then calibrated.

STANDARDS REQUIREMENTS:

- For scales of 7500g: 5 kg and 7.5kg
- For scales of 10000g: 5 kg and 10kg

V. TO RESPECT

1. Use the scale only as explained above in the technical details.
2. The unit should be leveled using the built-in spirit level.
3. Use the device on a stable and flat surface. Protect the device from shocks, humidity, radiations and magnetic fields.
4. If the screen does not show a stable result, make sure the support of the platform is not damaged.

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