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für den

## Pipettierhelfer witoped XP

### Einführung

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for the  
**witoped XP Pipet-Aid**

## Introduction

The witoped XP is suitable for use with all normal graduated and volumetric pipettes up to 100 ml. It makes pipetting not only safer but, particularly for difficult media, problem-free.

## Important accessories and spare parts

Filter (5 pcs)	Order no. 5380.006
Rubber insert (2 pcs)	Order no. 5380.010
Check valve with button assembly and button spring	Order no. 5380.007
Stand	Order no. 5380.003

## Important safeguards

When using electrical appliances, basic safety precautions should always be followed. Please read all instructions carefully before using unit.

### To reduce the risk of fire, electrical shock or injury to persons:

- To protect against the risk of electrical shock, do not put unit or charger in water or other liquid.
- Do not operate in the presence of explosives and/or flammable fumes or liquids.
- Do not operate charger with damaged cord or if damaged in any way.
- Use charger only in a standard electrical outlet.
- Do not attempt to use the charger with any other product. Do not attempt to charge this unit with any other Charger.
- Do not handle plug of charger with wet hands.
- This product includes a rechargeable NiCd 9V battery. Do not incinerate batteries as they will explode at high temperatures.

## Operating instructions

Insert a pipette into the nosepiece by twisting the pipette when pushing. Exercise extreme caution when using glass as breakage is an ever-present danger. Depressing the top button fills the pipette, depressing the bottom button empties it. The flow rate is proportional to the amount the button is depressed.

## Speed control

Each button has three speeds: "F" fast, "M" medium, "S" slow. The speed is selected by rotating the button on its axis until a detent is felt; the speed chosen is visible when the handle is held in the right hand. The "M" corresponds to the speed of previous Pipet-Aids. Fast is to be used for rapid filling of large pipettes or breaking up cell clumps. Slow speed should be used for small pipettes or where precise delivery is required. Please note that each button can be set independently so, for example, one could fill a large pipette on "F" and deliver small amounts on "S".

### Caution !

Speed control buttons must be positioned on "F", "M" or "S".  
Intermediate settings will cause leakage.

Do not change speed control buttons with fluid in the pipette. This will also cause the unit to leak.

## Battery

The XP is powered by a NiCd rechargeable battery (9V, 120 mAh) which must be disposed of in a proper manner after use. To get the maximum life from the battery, recharge it overnight as needed (recharging time about 15 h). The battery can be replaced by removing the screw marked (B) on diagram and sliding out the cover (78).

## Filter (0,8 µm, non-sterile)

The XP incorporates a TISSUE CULTURE NOSEPIECE, which features a unique filter element. The TC filter will cease to pass air in either direction when it becomes wet, alerting the user to possible contamination of the rubber insert. To remove the filter, slide the nosepiece release (74) away from the nosepiece and twist off the nosepiece.

## Components

Included in this brochure is a parts list and assembly drawing. Commonly used parts (e.g. rubber inserts, battery) are available from your dealer or witeg.

## Warranty

A quality product earns respect. This is why we provide a warranty period of 6 months for the witeped XP; this covers material and manufacturing errors.

We assume no responsibility for damage resulting from incorrect handling.

The instrument is carefully checked before leaving the factory. Nevertheless it is not possible to completely avoid faults occurring in use.

Checking returned instruments takes a lot of time and is expensive. This is why we require you to provide us with details of the faults. If we receive returned components without such details we will apply an additional processing charge of 25,- DM.

### To process your complaints we require the following details:

(please photocopy, fill in and return with the instrument)

Serial number: \_\_\_\_\_

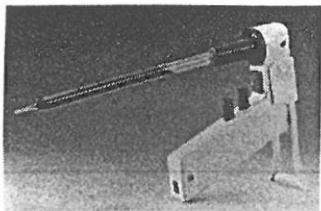
Purchase date /dealer: \_\_\_\_\_

Type of fault: \_\_\_\_\_

\_\_\_\_\_  
\_\_\_\_\_

Type of sample: \_\_\_\_\_


Sample temperature: \_\_\_\_\_

## Instrument stand

Included with your witeped XP is a convenient slide-on stand which allows you to set the Pipet-Aid down without removing or emptying the pipette. This stand will prevent liquid from running back into the nosepiece. When not needed, simply slide it off and store in a convenient place.

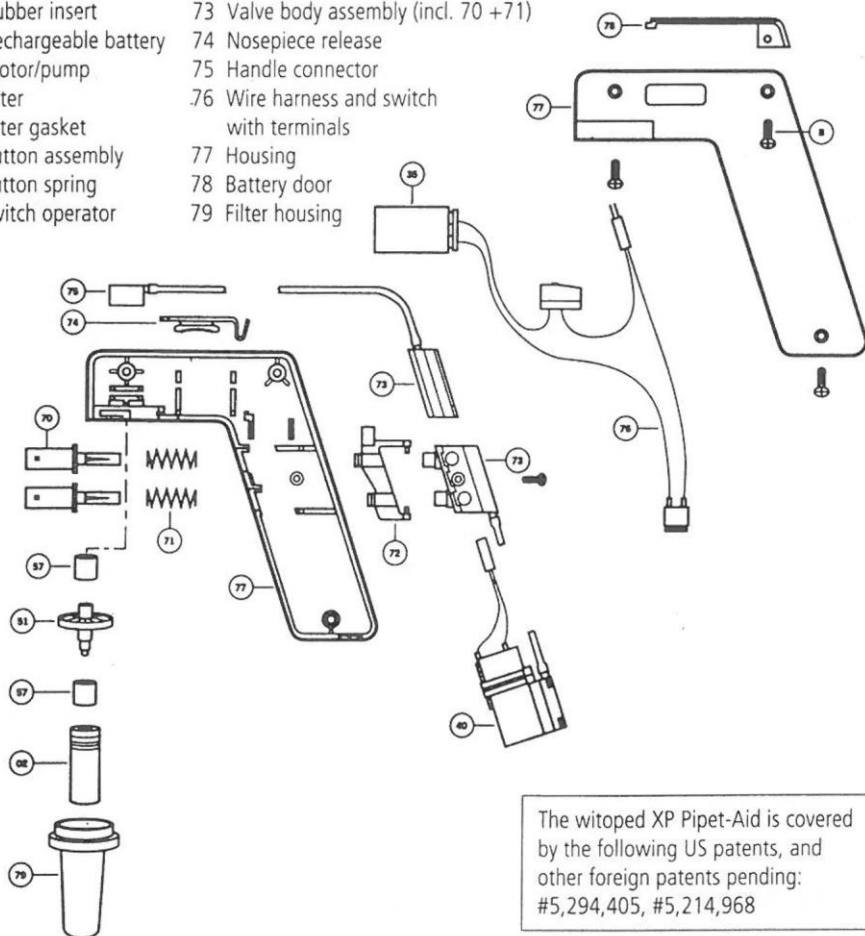- |                         |   |
|-------------------------|---|
| 02 Rubber insert        | 73 Valve body assembly (incl. 70 +71)     |
| 35 Rechargeable battery | 74 Nosepiece release                      |
| 40 Motor/pump           | 75 Handle connector                       |
| 51 Filter               | 76 Wire harness and switch with terminals |
| 57 Filter gasket        | 77 Housing                                |
| 70 Button assembly      | 78 Battery door                           |
| 71 Button spring        | 79 Filter housing                         |
| 72 Switch operator      |   |

The witeped XP Pipet-Aid is covered by the following US patents, and other foreign patents pending:  
#5,294,405, #5,214,968

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**Witeg Labortechnik GmbH**