ISO 9001  
CERTIFIED

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## **Instruction Manual**

*Please read before use!*

# **pipetus®-junior**

*manuelle Pipettierhilfe*

*Manual pipetting aid*

P990 30 01

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## *Contents*

	<i>Page</i>
1. Safety instructions	3
2. Operating principle	6
3. Pipetting	7
4. Maintenance	8
5. Available spare parts	8

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## 1. Safety Instructions

**Attention:** Before use please read the safety instructions and observe them carefully. Before use every member of staff must be familiar with these safety guidelines and they must be kept available at all times. This instruction manual cannot, however, cover every possible safety hazard. It is every user's responsibility to observe all health and safety precautions and to determine and observe restrictions before use of the unit.

- Observe all general safety practices and guidelines for laboratories, such as the use of protective clothing, safety glasses and gloves when handling hazardous chemicals.

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**Attention:** If the unit is opened, the warranty is invalidated and any resultant claims void.

- The filter supplied contains a hydrophobic 0,2 µm membrane for air filtration in either direction.
- Only have the unit professionally serviced.
- Never use undue force!

**Attention:** Only ever use the original accessories and replacement parts to avoid damaging the unit.

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## **2. Operating principle**

*pipetus®-junior* has no necessity of an electrical power supply, as a manually operated suction pump is built-in the handel. *pipetus®-junior* is suitable for all pipettes from 0,5 ml upwards. Pipetting is controlled by separate fingertip control buttons to fill (PUMP) and dispense (OUT). Pipette filling is controlled by repeated pumping of the button, with each pump about 2 ml of liquid are aspirated into the pipette. Dispensing is controlled by a control button (OUT). By pressing the OUT button the pipette will deliver by gravity. When using blow-out pipettes, the residual liquid in the pipette tip is ejected through unique pressure on the rubber ball (A). The speed of delivery can be regulated by pressing the OUT button.

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### **3. Pipetting**

Insert a pipette into the pipette holder. Fill the pipette to slightly above the required volume by pressing the pump button. Adjust to the exact volume by gently pressing the OUT button. To empty the pipette, press the OUT button. If using blow-out pipettes, the residual volume can be blown-out by pressing the rubber ball (A).

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#### **4. Maintenance**

When normally using the unit, it is maintenance free. Depending on the use, the filter (99 37 04 - set of 5 pcs) will require replacement. The complete nose piece assembly (990 23 00), except the disposable filter, can be sterilized by autoclaving (121°C).

#### **5. Available spare parts**

990 23 00	pipetholder junior complete
990 37 01	pipetholder case
990 37 04	round filter set with one-sided connection (5 filters)
990 32 01	silicone adapter
990 20 37	spacer